

AGTRON

Procedure:

- 1) To turn on machine, press the red button on the front panel, warm up for 30 minutes prior to first reading.
- 2) Pull drawer open and adjust calibration knob so that the needle is aligned with the calibration mark
(NOTE: Calibration mark ~48)
- 3) The Agtron and LED use the same sample dish and volume. Make sure the sample is smooth and with no bubbles.
- 4) Carefully place it into the drawer and close.
(NOTE: Make sure the tray is clean before placing the sample dish in the machine. This will avoid errors)
- 5) Allow a few seconds for the needle to settle then record